

# Purification of 6xHis AdK-oligo

Protein folding  
Protocol  
2024



# Protocol introduction

## Purification of 6xHis AdK-oligo using His SpinTrap™ (GE28-4013-53)

### Notes before starting the purification

- Modifications applied to the instructions of the manufacturer (SpinTrap™) are underlined and are recommended to improve the yield of the protein-oligo purifications.
- It is highly recommended to follow handling instructions of the manufacturer, including safety warnings and precautions.

### Buffer preparation

**Note:** Make sure that the imidazole solution used to prepare the following buffers has a pH ranging between 7 and 8.

- Prepare binding buffer: 40 mM Tris-HCl pH 7.5, 300 mM NaCl, 20 mM imidazole.
- Prepare elution buffer: 40 mM Tris-HCl pH 7.5, 300 mM NaCl, 500 mM imidazole.

### Column equilibration

- Following the manufacturer's instructions:
  1. Invert and shake the column repeatedly to resuspend the medium containing the His SpinTrap™ beads (resin)
  2. Loosen the top cap one-quarter of a turn and twist off the bottom closure.
  3. Place the column in a 2 ml microcentrifuge tube and centrifuge for 30 s at 70 to 100 x g.
  4. Discard the flow-through.
  5. Remove and discard the top cap.
- Add 600 ul of binding buffer to the resin in the column and mix by pipetting with a cut tip until the resin is homogeneously resuspended.
- Centrifuge for 30 s at 70 to 100 x g, discard flow-through, and re-use the 2 ml tube for the next step.

---

## Binding of 6xHis AdK-oligo to the resin

- Upon completion of the Step 1 “Purification of protein-oligo chimera” of Protein labeling (cysteine or ybbR), add 80  $\mu$ l of binding buffer to the 80  $\mu$ l of 6xHis AdK-oligo (Cys) solution or 115  $\mu$ l of binding buffer to the 45  $\mu$ l of 6xHis AdK-oligo (ybbR). **Note:** all the following steps are identical for 6xHis AdK-oligo (Cys) and 6xHis AdK-oligo (ybbR).
- Add the 6xHis Adk-oligo solution diluted with binding buffer to the resin and mix by pipetting with a cut tip until the resin is homogenously resuspended.
- Incubate for 3 minutes. Then resuspend the resin again homogenously.
- Repeat the previous step for additional two times to obtain a total incubation time of AdK-oligo with resin of  $\sim$ 9-10 min. **Note:** if a shorter incubation time is preferred for the protein of interest, go to the next step after the first 3 minutes incubation with the resin.
- Centrifuge for 30 s at 70 to 100 x g, discard flow-through, and re-use the 2 ml tube for the next step.

## Clean up from the excess of unreacted oligos

- Add 600  $\mu$ l of binding buffer to the resin and resuspend the resin homogenously.
- Centrifuge for 30 s at 70 to 100 x g, discard flow-through, and re-use the 2 ml tube for the next step.
- Repeat the previous two steps for three additional times to ensure a complete clean up from the excess of unreacted oligos.

## Elution

- Place the column in a new 2 ml tube.
- Add 200  $\mu$ l elution buffer and resuspend the resin homogenously.
- Incubate for 3-5 min.
- Centrifuge for 30 s at 70 to 100  $\times$  g.
- Using the same 2 ml tube, add 200  $\mu$ l elution buffer and resuspend the resin homogenously.
- Incubate for 3-5 min.
- Centrifuge for 30 s at 70 to 100  $\times$  g and collect the purified 6xHis AdK-oligo protein.

You are now ready to continue with the buffer exchange from the first bullet point in Step 2 Purification of protein-oligo chimera.

---

[info@lumicks.com](mailto:info@lumicks.com)  
[www.lumicks.com](http://www.lumicks.com)

Or find us on:



### LUMICKS HQ

Paalbergweg 3  
1105 AG Amsterdam, The Netherlands  
+31 (0)20 220 0817



### LUMICKS USA

800 South Street, Suite 100,  
Waltham, MA, 02453, USA  
+1 781 366 0380



### LUMICKS Asia

Room 545, Block A, Langentbldg Center  
No.20 East Middle 3rd Ring Road  
Chaoyang District, Beijing, 100022 China  
+86 (0) 10 5878 3028

All content and images used in this document are owned or licensed by LUMICKS B.V. Unauthorized use is prohibited. Any information provided herein by LUMICKS is made available "as is" and [you] understand and agree that such information is made available without any representation or warranty, express or implied, including any implied warranty of merchantability, satisfactory quality or fitness for any particular purpose or any warranty that the use of such information will not infringe or violate any patent or other proprietary rights of any third party. For the latest product information please consult us directly.

C-Trap®, m-Trap®, AFS®, u-Flux™, Bluelake™, z-Movi®, LUMICKS and the LUMICKS logo are registered trademarks of LUMICKS B.V.

© LUMICKS B.V. Amsterdam, The Netherlands.

# LUMICKS